

# UPB-150-ART

<https://www.gigahertz-optik.com/en-us/product/upb-150-art/>


**Product tags:**



## Description

### Photometry to determine material properties

Material properties is a general term used to refer to the reflection, transmission and absorption properties of a material. Photometry hereby refers to a measurement process through which these properties are determined using a photometer.

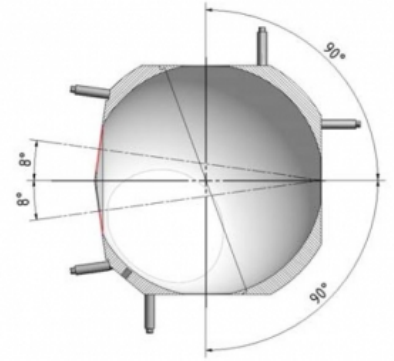


*Positions of the five measurement ports of the UPB-150-ART enable its universal application.*

### Integrating sphere photometer

The integrating sphere is very common in both broadband and narrow-band photometric applications for diffuse sample illumination and determination of the reflected and transmitted light proportions.

### UPB-150-ART precision integrating sphere

With the 150mm diameter, the UPB-150-ART provides a good compromise in terms of precision of the sphere function and level of the measurement signal. Its manufacture from an aluminum block ensures a very precise alignment of the measurement ports to each other. Another characteristic property are the thin circumferential knife-edges of the measurement ports that enable a high acceptance angle for diffuse light.

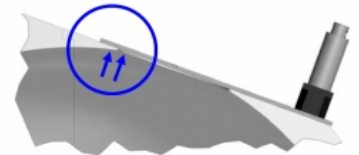


*The front side of the port plugs, port reducers and sphere surface are at the same level and thus ensure an almost ideal reflection surface.*

### Multifunctional integrating sphere

The UPB-150-ART is ideal for multiple functions through its five knife-edged measurement ports that can be used either while open or closed. Possible measurement setups include:

- 0°/d Reflection
- 8°/d reflection with and without glamour
- 0°/d directed and diffuse transmission
- Measurements with correction of sample absorption effects
- ~~Two-way measurements~~

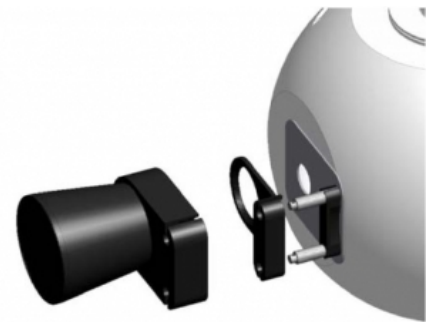

*Light trap for absorption of the transmitted light of transparent samples in 8°/d reflectance measurement setup. Combined with port reducer.*

### Accessories for individual sphere assemblies

Port plugs and port reducers that can easily be fixed and removed using the Slide & Fix holder are offered for either closure or reduction of ports. An important feature of the UPB-150-ART is that the front side of the plugs and reducers projects into the port and is therefore on the same level as the sphere surface.

### Universal sample holder

The sample holder is fixed onto the posts of the Slide & Fix holder and ensures secure mounting of thin and thick samples. An important feature of the Slide & Fix system is that the sample holder can be combined with port reducers and a light trap.



*BTS2048-XX-F spectrometer used as detector*

---

## Light trap with Slide & Fix holder

The light trap has a 30mm input port and ensures minimal back reflection over the entire usable spectral range of the integrating sphere through its internal design and black absorbing coating. The Slide & Fix holder can be used to combine the light trap with port reducers and the sample holder.

---

## Ex-changeable holders for detectors and light guides

Ex-changeable adapters are offered for mounting of detectors and light guides. Each has a baffle that can be aligned to one of the five measurement ports accordingly depending on the measurement application. The UMPA port adapter of our modular integrating spheres can hereby be used. It is also possible to connect one of our fiber-coupled spectrometers [BTS2048-VL-TEC-F](#) or [BTS2048-UV-F](#).



*UPB-150-ART-Z01, port plug with flat surface*



*UPB-150-ART-Z02-3, 3mm port reducer*

---

## Reflection coating made of barium sulfate

The UPB-150-ART is coated with Gigahertz-Optik's self-manufactured diffusely reflective barium sulfate (ODP97) coating. The multi-layer coating ensures good diffuse reflective properties over the entire usable spectral range.



*UPB-150-ART-Z02-15, 15mm port reducer*



*UPB-150-ART-Z02-35, 25mm port reducer*



*UPB-150-ART-Z03, sample holder*



*UPB-150-ART-Z04, light trap*



*UPB-150-ART-Z05, 0.5" detector adapter with baffle*



*UPB-150-ART-Z06, 1.0" detector adapter with baffle*



*UPB-150-ART-Z07, 2.0" detector adapter with baffle*



*UPB-150-ART-Z09, port plug with spherical surface*



*UPB-150-ART-Z10, port reducer with spherical surface. Free diameter if specified*

## Specifications

### General

Short description	Five port integrating sphere for universal use in transmission and reflection measurement application in the 300 nm to 2500 nm spectrum.
Main features	Barium sulfate coating. Machined out of solid material. Therefore very precise spherical shape, position of the application ports and finish of the knife edges of the application ports. Accessory for the application ports with Slide&Fix mounting System. Detector port with baffle.
Measurement range	Diameter: 150 mm (6 inch) Coating: Barium sulfate. Spectral range 300 nm to 2500 nm Application ports: five, 30 mm diameter with knife edge, post for Slide&Fix accessory Detectorport: exchange able Adapters with baffle
Typical applications	Spectrophotometer. Other transmission and reflection measurement set-ups.
Calibration	-
<b>Product</b>	
Sphere diameter	150 mm


Coating	ODP97 (Barium Sulfate)
Measurement Port Diameter	30 mm, with knife edge design
Knife edge thickness	0.5 mm
Number of measurement ports	Five
Angle tolerance of port setup	$\pm 0.1^\circ$
Accessories	Two detachable posts for the Slide & Fix system per port
Detector port	30 mm with circumferential frame ( $\varnothing = 58$ mm) with lateral V-notch
Mounting	Flat surface with four M5 threaded bores on a 50 mm pitch circle
<b>Miscellaneous</b>	
Weight	approx. 1.6 kg

## Downloads

Type	Description	File-Type	Download
Drawing	UPB-150-ART	pdf	<a href="https://www.gigahertz-optik.com/assets/Uploads/102257_UPB-150-ART.pdf">https://www.gigahertz-optik.com/assets/Uploads/102257_UPB-150-ART.pdf</a>

## Configurable with

Product Name	Product Image	Description	Go to product
UFC-11		Fiber adapter for use with UP and UM type integrating spheres. Features: SMA, FC, SC and ST connectors.	<a href="https://www.gigahertz-optik.com/en-us/product/ufc-11/">https://www.gigahertz-optik.com/en-us/product/ufc-11/</a>
VL-1101		Photometric detector head with VL-11 mount. Features: modular detector for use with integrating spheres, front lenses etc. For use with optometers and signal amplifiers	<a href="https://www.gigahertz-optik.com/en-us/product/vl-1101/">https://www.gigahertz-optik.com/en-us/product/vl-1101/</a>
BN-Rxx-D2		Calibration Standards for Diffuse Reflection	<a href="https://www.gigahertz-optik.com/en-us/product/bn-r98-d2/">https://www.gigahertz-optik.com/en-us/product/bn-r98-d2/</a>
ODMP		Synthetic Diffuse Reflecting Material in Different Gray Scales	<a href="https://www.gigahertz-optik.com/en-us/product/odmp/">https://www.gigahertz-optik.com/en-us/product/odmp/</a>

Product Name	Product Image	Description	Go to product
CLS-21		Collimator Light Source for Integrating Spheres	<a href="https://www.gigahertz-optik.com/en-us/product/cls-21/">https://www.gigahertz-optik.com/en-us/product/cls-21/</a>

## Purchasing information

Article-Nr	Modell	Description
<b>Product</b>		
15297566	UPB-150-ART	Integrating sphere
<b>Accessories</b>		
15297567	UPB-150-ART-Z01	Port plug
15297568	UPB-150-ART-Z02	Port reducer. xx aperture diameter set in 0.1mm steps.xx =
15297496	UPB-150-ART-Z02-3	Port reducer 3mm
15297497	UPB-150-ART-Z02-15	Port reducer 15mm
15297498	UPB-150-ART-Z02-25	Port reducer 25mm
15297569	UPB-150-ART-Z03	Sample holder
15297570	UPB-150-ART-Z04	Light trap
15297571	UPB-150-ART-Z05	UMPF-0.5 Detector adapter
15297572	UPB-150-ART-Z06	UMPF-1.0 Detector adapter
15297573	UPB-150-ART-Z07	UMPF-2.0 Detector adapter
15298198	UPB-150-ART-Z09	Port plug
15298199	UPB-150-ART-Z10	Port reducer. xx aperture diameter set in 0.1mm steps.xx =

## Contact, Calibration, Service & Support

We are known worldwide for excellent technical consulting and after sales support. Contact us to find together the best solution for you. Our services:

- Technical Consulting & Sales
- After-Sales Support
- Calibrations & Re-Calibrations ([ISO/IEC 17025 Calibration Services](#), [factory calibration](#), [Calibration of Third-Party Products](#))
- Repairs & Updates
- OEM & Feasibility Consulting of Customized Solutions

[Send us your inquiry](#) or contact us by phone or e-mail. We would welcome your feedback too or review us on [Google](#).

### Gigahertz Optik GmbH (Headquarter)

Tel.: +49 (0)8193-93700-0  
Fax: +49 (0)8193-93700-50  
[info@gigahertz-optik.de](mailto:info@gigahertz-optik.de)

An der Kaelberweide 12  
82299 Tuerkenfeld, Germany

### Gigahertz-Optik, Inc. (US office)

Phone: +1-978-462-1818  
[info-us@gigahertz-optik.com](mailto:info-us@gigahertz-optik.com)

Boston North Technology Park  
Bldg B - Ste 205  
Amesbury, MA 01913 USA